

STEP 1

THE P-20 MEASURES 2-20 LIL, THE P-200 MEASURES 20-200 LIL

THE TOP SQUARE SHOWS 10'S IN THE P-20 AND 100'S IN THE P-200

STEP 2

THE BOTTOM SQUARE SHOWS .1'S (10THS) IN THE P-20 AND 1'S IN THE P-200

WHICH PIPET DO WE USE FOR 10 MICROLITERS?

SELECT THE CORRECT PIPET FOR THE AMOUNT YOU NEED

DIAL THE AMOUNT YOU NEED ON THE PIPET

STEP 3

PUT A TIP ON THE PIPET

CLOSE THE TIP BOX QUICKLY!

THAT'S RIGHT, LET'S KEEP THEM STERILE!

STEP 4

YEAH! AND KEEP ALL THE LIQUID IN THE TIP, NOT IN THE PIPET.

KEEP THE PIPET VERTICAL

STEP 5

I NEED TO HOLD THE PIPET AND THE TUBE MYSELF!

PLACE PIPET AND SOURCE TUBE AT EYE LEVEL

STEP 6

WAIT! DON'T PRESS IT WITH THE TIP INSIDE THE TUBE!

PRESS THE PULLER TO THE 1ST STOP

STEP 7

STEP 8

STEP 9

STEP 10

STEP 11

IF I PUT TOO MUCH OF THE TIP IN, MORE LIQUID WILL BE WASTED ON ITS SURFACE.

FOLLOW THE SURFACE OF THE LIQUID DOWN WITH THE TIP.

PUT THE TIP JUST BELOW THE SURFACE OF THE LIQUID

RELEASE THE PULLER SLOWLY

HAND ME THE CLEAN TUBE AND TAKE THIS ONE, PLEASE! I STILL HAVE TO HOLD THE TUBE MYSELF.

PUT TIP IN TARGET TUBE

IF I PRESS SLOWLY, I MAY NOT NEED TO PRESS TO THE SECOND STOP

DON'T LET GO OF THE PULLER UNTIL YOU TAKE THE TIP OUT OF THE TUBE!

PRESS TO 2ND STOP AND REMOVE TIP FROM TUBE

WE DID IT! GOOD JOB!

RELEASE PULLER SLOWLY AND THROW TIP IN WASTE.